

# Blades Biological Ltd

COWDEN . EDENBRIDGE . KENT . UK . TN8 7DX

Telephone: (01342) 850242 . Fax: (01342) 850924

Email: [info@blades-bio.co.uk](mailto:info@blades-bio.co.uk) Online: <http://www.blades-bio.co.uk/>

## **Alga-gro Concentrated Medium**

Alga-gro Concentrate is a sterile enrichment for both seawater and freshwater, producing an almost universal algal medium.

### **Freshwater Medium**

For each litre of medium required, add 20ml tube to 1 litre of bottled mineral water.

If sterile medium is required, transfer into containers and autoclave at 15lb pressure for 15 minutes.

### **Seawater Medium**

Seawater should be sterilized before use either by boiling or autoclaving at 7lb pressure at 100c for 3 minutes.

When cold, mix 20ml tube of alga-gro to each litre of seawater required.